



Explorer[®] Balance Display Tower Mount Kit

Installation Instructions

1. INTRODUCTION

Please read this manual completely before installation and operation.

CAUTION:



READ ALL SAFETY WARNINGS BEFORE INSTALLING, MAKING CONNECTIONS, OR SERVICING. FAILURE TO COMPLY WITH THESE WARNINGS COULD RESULT IN PERSONAL INJURY AND/OR PROPERTY DAMAGE. RETAIN ALL INSTRUCTIONS FOR FUTURE REFERENCE.

1.1 Kit Contents



Five M5 Screws

Hex Wrench

Tower with Top
and Foot

2. INSTALLATION

2.1 Safety Precautions

For safe and dependable operation of this equipment, please comply with all safety precautions mentioned in this manual.

Disconnect equipment from mains power supply before installation.

2.2 Installation procedure


Follow the steps below. Refer to the Figures when performing these steps.

1



Make sure Balance is unplugged (turned OFF).

2

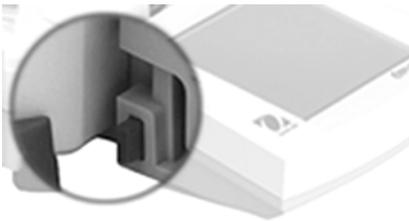

Detach the Display:
press buttons on both sides to disconnect. Then pull Display away from base.

3




Open the Display:
press buttons (enlarged in photo),
on both sides, then lift Display until
it opens.

4



Open the Display to
access and release the
Display Communication Cable.

5

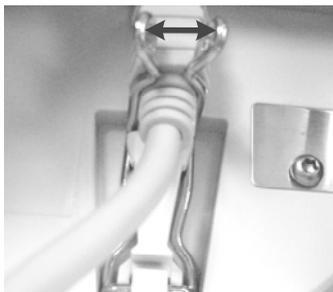

Release the clip that secures the Display Communication Cable: press out on both sides, as shown.


Unplug the Display Communication Cable.

6



Un-spool the Display Communication Cable from the base.

7

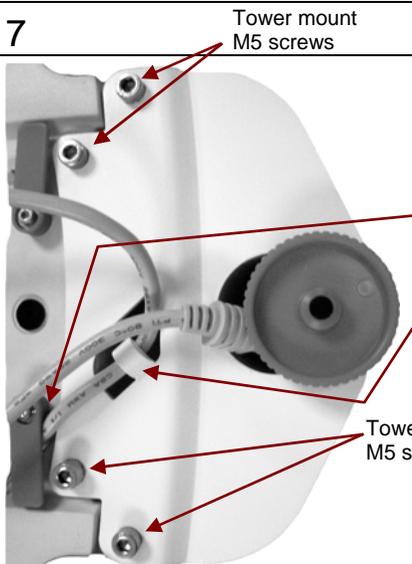


Attach Tower to the Base using the four M5 screws and wrench (supplied).


Secure the cable as shown:

1. First, under the cable guide,
2. Then under the groove in the tower base.

Insert the cable plug in the tower base and feed it through the tower.

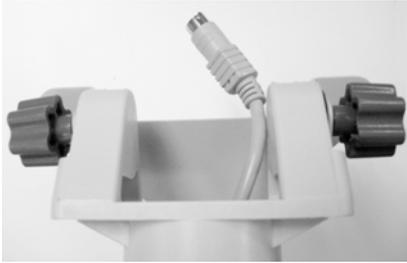8


Pull the Display Communication Cable through the top of the tower.

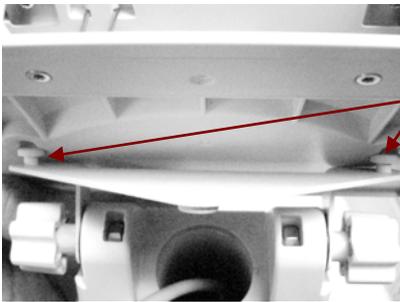9

Position the Display above the Tower Platform.

Insert the posts in the receptacles in the Display Base.

Slide the Display forward as shown, to seat it in the Tower platform.

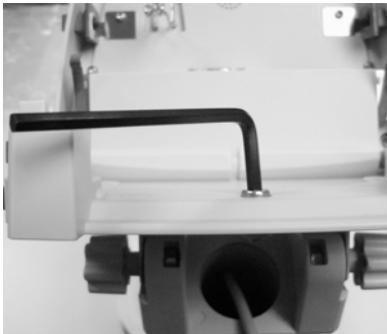10

Insert one M5 screw in the Display Base to secure it to the Tower Platform.

Use the supplied Hex Wrench to tighten the screw.

11



Insert the cable plug with the security clip open, then press the security clip into position to secure the cable.

Route the cable through a groove in either side of the Display, or from the rear of the Display, as preferred.

Pull excess cable down into the Tower for safety. (Best practice: remove slack and re-wrap cable around cable guides on the bottom of the balance.)

12



Re-connect power to the Base. Use the Level Assist feature to re-level the Balance. (Keep the Tower foot raised during this procedure.)

After leveling the Balance, adjust the Tower foot so it supports the Tower and is level with the feet of the Base.

13



To adjust the Display viewing angle:

- Press in the two Display Adjust buttons and move the Display.

Or...

- Loosen the knob that secures the Tower top and move it.



Ohaus Corporation
7 Campus Drive
Suite 310
Parsippany, NJ 07054, USA
Tel: (973) 377-9000
Fax: (973) 593-0359

With offices worldwide

www.ohaus.com



P/N 83021181 © 2011 Ohaus Corporation, all rights reserved.

Printed in China